

SOC Medium

Cat No: 40021, 40022

Storage: Room Temperature

Compositions:

2% Peptone, 0.5% Yeast extract, 10mM NaCl, 2.5mM KCl,
10mM MgCl₂ , 20mM MgSO₄, 20mM glucose

Operation procedures:

It is best to use all of the Medium once it is opened.

If the Medium has to be used multiple times after it is opened,
sterilizing operation is needed.

Wash both hands and clean the working surface by 70% ethanol.

Do not cough, sneeze, or breath in the media. Otherwise the
Medium will be contaminated.

Hold breath when opening the Medium and quickly close
it after transferring the Medium out.

Avoid opening the Medium in places such as bacterial culture
rooms.

Use a biological hood if contamination persists.

Warranty:

We guarantee the sterilization and composition of
our SOC Medium. If the product fails to meet its
specs, we will promptly replace the product. All
other warranties are expressly disclaimed.