

Sterilized LB Lennox Broth

Cat No: 40011, 40012

Storage: Room Temperature

Compositions:

1% Peptone, 0.5% Yeast extract, and 0.5% NaCl

Operation procedures:

1. It is best to use all of the medium once it is opened.
2. If the medium has to be used multiple times after it is opened, sterilizing operation is needed.
3. Wash both hands and clean the working surface by 70% ethanol.
4. Do not cough, sneeze, or breath in the media. Otherwise the medium will be contaminated.
5. Hold your breath when opening the medium and quickly close it after transferring the medium out.
6. Avoid opening the medium in places such as bacterial culture rooms.
7. If you still cannot avoid contamination, you need a biological hood.